

ScanR Switch On/Off Procedures

1. Check that the computer is ON.
2. Switch on ScanR (Wall plug labelled ScanR on the lefthand side of the microscope frame)
3. Switch on camera (Wall plug labelled Camera on the left-hand side of the microscope frame)
4. Switch on microscope controller (on the left-hand side of the microscope frame on the bench. Switch is on the back panel)
5. Log into your university account
6. Run ScanR Acquisition

1. Close ScanR Acquisition
2. Log off
3. Switch off camera (Wall plug labelled Camera on the left hand side of the microscope frame)
4. Switch off ScanR (Wall plug labelled ScanR on the left-hand side of the microscope frame)
5. Remove your slides and clean the stage

Need help? Contact Yan Gu: 07981875417, office: 2c37, ext:6997