

ScanR_Robot Switch On/Off Procedures

1. Switch on microscope (Wall plug labelled Microscope on right-hand of the chamber)
2. Switch on camera (Wall plug labelled Camera on right-hand of the chamber)
3. Switch on robot loader (Only when needed! Wall plug labelled Loader on left-hand of the chamber)
4. Switch on microscope controller (Sitting on bench, on right-hand side of the chamber labelled Controller. Switch is at the back panel)
5. Log in
6. Run scanR Acquisition software (When the controller finishes initialization)

1. Close scanR Acquisition software
2. Log off
3. Switch off camera (Wall plug labelled Camera on right-hand of the chamber)
4. Switch off microscope (Wall plug labelled Microscope on right-hand of the chamber)
5. Switch off robot loader (Only when the loader is switched on! Wall plug labelled Loader on left-hand of the chamber)
6. Shut all chamber doors