

# STED Switch On/Off Procedures

1. Make sure the workstation is on
2. Switch on STED control (wall plug labelled STED behind screen)
3. Switch on microscope touch pad (back side of touch pad)
4. Log into university account
5. Wait until initialization of touch pad finishes
6. Start Inspector Lightbox software

You may need to turn on the microscope 3hrs in advance to stabilize the microscope.

1. Close Impector software
2. Log off university account
3. Switch off STED control (wall plug labelled STED behind screen)
4. Take off the sample
5. Clean the objective (use lens tissue only)

**Need help?** Contact Yan Gu: 07981875417, office: JMS 2c37