

AxioVert Switch On/Off Procedures

1. Check that the computer is ON.
2. Switch on microscope (Wall plug labelled Stripe 3)
3. Switch on camera (camera controller labelled Camera on shelf)
4. Switch on fluorescence lamp (HXP120V lamp box on shelf labelled Lamp)
5. Log into your university account
6. Run Micro-manager

1. Close Micro-manager
2. Log off
3. Switch off fluorescence lamp (HXP120V lamp box on shelf labelled Lamp)
4. Switch off camera (camera controller labelled Camera on shelf)
5. Switch off microscope (Wall plug labelled Stripe 3)
6. Remove your slides and clean the objective with lens tissue

Need help? Contact Yan Gu: 07981875417, office: 2c37, ext:6997